

Technical Note

Constructing an interface cable between the Kistler Quattro Jump force plate and BMS XP6 Interface Unit.

Most Recent Revision Date: 16 June 2004

1. The cable consists of:
 - a. 2-3 meters of 2 core shielded cable
 - b. 1 x male DB15 plug and backshell
 - c. 1 x 8 pin RJ45 socket



2. Wire in one end of the cable to the cable to the RJ45 with one wire to terminal 1 and the other to terminal 6.



3. Solder the DB15 on to the other end of the cable. The connections are:



